

## Normal Donkey serum blocking buffer

#Cat: NB-23-00183-1

Size: 100ml

#Cat: NB-23-00183-2

Size: 18ml

### Intended Use:

Normal Donkey Serum Blocking Buffer is used in immunohistochemical staining procedures to block non-specific staining when the host species of the secondary antibody is Donkey.

### Formulation:

Solution contains 0.05% NaN<sub>3</sub>.

### Recommended Protocol

1. Apply 100uL or enough volume of Normal Donkey Serum Blocking Buffer on tissue before adding primary antibody and incubate for 10 minutes at room temperature.
2. Blot off Normal Donkey Serum Blocking Buffer Solution (Do not rinse) and add primary antibody.

### Storage

Store at 2-8°C. Return the reagent back to 2-8°C after each use.

### Precaution:

Wear gloves and take other necessary laboratory safety procedure

*For reference only*

*For Research Use Only. Not for Diagnostic or Therapeutic Use.*